

Precision at your fingertips



STEPPER™

REPEATER PIPETTE FOR ECOSTEP™ SYRINGES



Stepper™ 416 repeater pipette

Latest generation of a compact and reliable instrument intended for serial dispensing within 10 to 5000 µL. Associated with three Ecostep™ syringes only, it provides a broad choice of 53 different volumes, and up to 73 doses per filling.

Stepper™ has a unique trigger action mechanism eliminating thumb fatigue. It enables a precise activation movement and is best suited for repeat dispensing, aliquoting, tubes and plate filling. Two-year warranty.

Main values

- 4-finger friendly trigger activation
- Setting knobs showing volumes and number of aliquots
- Only three positive displacement syringes
- Colour coded setting knobs and syringes
- Self-locking mechanism
- Regular and *bioproof*™ Ecostep™ syringes
- Broad chemical compatibility



Instant volume reading ①

Select appropriate volume by simple rotation. Volume and corresponding number of aliquots are clearly visible.

Syringe filling ②

Pull trigger to aspirate liquid for full or partial syringe filling.

Ready to dispense ③

Instrument fits any hand in a very comfortable way.

Finger activation with no thumb fatigue ④

Ergonomic, effortless distribution movements using four fingers.

Release for next dose ⑤

Easy trigger action to monitor fast or slow dispensing.

Self-locking mechanism ⑥

Stop function prevents false delivery if remaining liquid is insufficient.